

# Perform Sterile Alcohol EP

## STERIL, ALKOHOLBASERET OPLØSNING

- Hurtig, effektiv og afsætter ingen rester
- Steril (gamma steriliseret)  
microbial filtreret (0,2 µm)
- Renrum grade A/B, ISO 5/6
- Klar til brug i handy og steril 500 ml sprayflaske
- Farve og parfumefri
- Effektivitetstest efter europæisk standard

## MIKROBIOLOGISK EFFEKTIVITET

- Bactericidal EN1276, EN13697  
Kontaktid: 1 min.
- Fungicidal EN1650, EN13697  
Kontaktid: 2 min.
- Verucidal  
Kontaktid: 30 sek.



VARENUMMER	TYPE	PAKNING
121802	500 ml sprayflaske	10 x 500 ml/karton



Sterile, alcohol-based ready to use solution for disinfecting surfaces in cleanrooms.

## perform® sterile alcohol EP

### Our Plus

- fast, effective and residue free
- sterile ( $\gamma$ -irradiated) and microbial filtrated (0.2  $\mu\text{m}$ )
- suitable for cleanrooms grade A/B, ISO class 5/6
- ready to use in a handy, sterile 500 ml spray system
- colour and perfume free
- efficacy tested according to European norms (EN)

### Application areas

Disinfection of surfaces and equipment/devices in cleanrooms. Especially for Grade A and B clean areas, according to:

- EU Guide to Good Manufacturing Practice for medicinal products Annex 1: Manufacture of sterile medicinal products, March 2009.
- FDA Draft Guidance for Industry: Sterile Drug Products produced by aseptic processing- current Good Manufacturing practice, U.S. Department of Health and Human Services; Food and Drug Administration, September 2004.

### Instructions for use

Apply perform® sterile alcohol EP undiluted to objects and surfaces and allow to take effect. Ensure complete wetting. The product dries off quickly. perform® sterile alcohol EP can be used for glove disinfection. Maximum application rate 50 ml/sq.m. Do not use to treat surfaces that are particularly sensitive to alcohol, such as acrylic glass.

Efficacy	Concentration	Contact time
bactericidal EN1276, EN13697	ready-to-use	1 min.
fungicidal EN1650, EN13697	ready-to-use	2 min.
virucidal against enveloped viruses in accordance with DVV (German Registered Association for Combating Viral Diseases)//RKI Guideline	ready-to-use	30 sec.



# perform<sup>®</sup> sterile alcohol EP

## Product data

Composition:

100 g solution contains the following active substances: 25 g Ethanol (94 %), 35 g Propan-1-ol.

## Chemical-physical data

Color	colourless
Density	ca. 0,89 g/cm <sup>3</sup> / 20 °C
Flash point	27 °C / Method : DIN 51755 Part 1
Form	liquid
pH	ca. 7 / 20 °C

## Special advice

Use disinfectants safely. Always read the label and product information before use.

The sterile 500 ml spray system offers the following advantages:

- convenient size for storage
- ergonomically shaped spray pump for ease of use
- remains sterile after opening

The bottles are sterile (with irradiation indicator) and doublewrapped individually. perform<sup>®</sup> sterile alcohol EP comprises alcohol of pharmaceutical quality in accordance with the Ph. Eur.

## Information for order

Item	Delivery form	Item no.
perform <sup>®</sup> sterile alcohol EP 500 ml bottle	10/Carton	on request

These products are not available in every country. For more information please contact our local subsidiary or distributor.

## Application aids

Application aids	Item no.
perform <sup>®</sup> spray bottle holder	693540
perform <sup>®</sup> spray bottle holder- plate	693541



Schülke & Mayr GmbH holds a Manufacturer's Authorisation according to sect 13 para 1 German Drug Law and Certificates of GMP Compliance for medicinal products.



A company of the Air Liquide-Group

## Related Products

- perform<sup>®</sup> sterile dry wipes
- perform<sup>®</sup> sterile mix dry wipes
- perform<sup>®</sup> sterile premium dry wipes

## Environmental information

schülke manufactures products economically and with advanced, safe and environmentally friendly production processes while at the same time maintaining out high quality standards.

## Expert opinion and information

Please visit our website for an overview of all available literature/reports on the product: [www.schuelke.com](http://www.schuelke.com).

For individual questions:

Customer Sales Service

Phone: +49 40 52100-666

E-Mail: [info@schuelke.com](mailto:info@schuelke.com)

schülke Headquarters  
Schülke & Mayr GmbH  
Robert-Koch-Str. 2  
22851 Norderstedt, Germany  
Phone +49 40- 52100- 0  
Fax +49 40- 52100- 318  
[www.schuelke.com](http://www.schuelke.com)  
[info@schuelke.com](mailto:info@schuelke.com)